

Comparison of Static and Dynamic Chemisorption Methods

1. Background

In previous issues of *AMI Notes*, we discussed the application of pulse chemisorption and temperature-programmed desorption (TPD) for catalyst characterization. A frequently asked question is how results from these dynamic methods compare with those obtained from static volumetric chemisorption. This note addresses that question by comparing data collected for various catalysts using both static and dynamic approaches.

2. Pulse Chemisorption

2.1 Pulsed Experiment

This technique is particularly well-suited for catalyst-adsorbate systems with relatively fast adsorption kinetics, meaning the adsorption process is not activated.

For this study, we selected two supported Pt catalysts with low metal loading:

- ✓ ASTM standard catalyst containing 0.5 wt% Pt on alumina
- ✓ In-house catalyst with 0.3 wt% Pt on alumina

Carbon monoxide (CO) was chosen as the adsorbate because CO adsorption is well-established as a rapid process, making CO the ideal choice for pulse chemisorption experiments. While hydrogen, CO, and occasionally oxygen are commonly used adsorbates for static (volumetric) chemisorption of Pt materials, hydrogen adsorption on Pt exhibits slow kinetics under dynamic conditions, making it a less suitable choice. ⁽¹⁻³⁾

2.2 Results

Table 1 presents a comparison of CO uptake measurements for the two Pt catalysts using both volumetric and pulse techniques. The results show excellent agreement between the methods. Notably, the pulse method yielded slightly lower uptake values, likely because the volumetric method also captures weakly held or "reversible" CO that the pulse technique does not detect.

Catalyst	Pt Loading (wt%)	Method	CO Uptake (μmol/g cat.)
ASTM Standard Pt/Al ₂ O ₃	0.5	Volumetric Chemisorption	10.2
ASTM Standard Pt/Al ₂ O ₃	0.5	Pulse Chemisorption	8.8
In-house Pt/Al ₂ O ₃	0.3	Volumetric Chemisorption	10.0
In-house Pt/Al ₂ O ₃	0.3	Pulse Chemisorption	9.1

Table 1: Comparison of volumetric and pulse CO chemisorption data on Pt/Al₂O₃

One significant advantage of the pulse chemisorption method over the volumetric approach is the time required to complete the analysis. After the necessary pretreatment, which is similar for both techniques, a full pulse chemisorption experiment typically takes less than 30 minutes, including system calibration. In comparison, a standard five-point volumetric measurement can easily require six hours or more.

A second advantage is the ease with which the sensitivity of the pulse method can be enhanced. This can be achieved by using smaller pulse loops or by diluting the adsorbate in an appropriate inert carrier gas. For example, a 10% CO in helium mixture can replace pure CO. While this may slightly increase the duration of the analysis, it significantly improves reproducibility.

3. Temperature-Programmed Desorption (TPD)

Temperature-programmed desorption (TPD) is especially useful when catalyst–adsorbate kinetics are not favorable for pulse chemisorption measurements. Cobalt metal catalysts are a typical example. For supported cobalt catalysts, hydrogen chemisorption proceeds slowly at room temperature but can be accelerated by raising the sample temperature to approximately 100°C.⁴ Carbon monoxide also chemisorbs slowly at room temperature and poses an additional challenge: at elevated temperatures, CO can disproportionate via the Boudouard reaction (Eqn. 1):



3.1 Experiment

In this study, the hydrogen uptake of two cobalt catalysts was measured using both volumetric chemisorption and TPD:

- ✓ 20 wt% Co/Al₂O₃
- ✓ 0.5 wt% Ru/20 wt% Co/Al₂O₃

The second catalyst contained added ruthenium as a promoter to aid reduction and improve dispersion.⁵

H₂ was chemisorbed onto the samples at 50 °C for 30 minutes, and then TPD was performed up to 400 °C. Static H₂ uptake was measured by a five-point volumetric method.

3.2 Results

Figure 1 presents the H₂ TPD profiles of these catalysts. Both exhibited broad desorption profiles, with the Ru-containing catalyst showing a larger desorption signal, as expected. The key parameters for obtaining reliable TPD results were an adsorption temperature of 50 °C and an adsorption time of 30 minutes.

Table 2 summarizes the H₂ uptake measured by both TPD and a five-point volumetric method. Once again, excellent agreement was observed between the two techniques. As with pulse chemisorption, TPD offers significant time savings compared to traditional volumetric methods. Furthermore, TPD

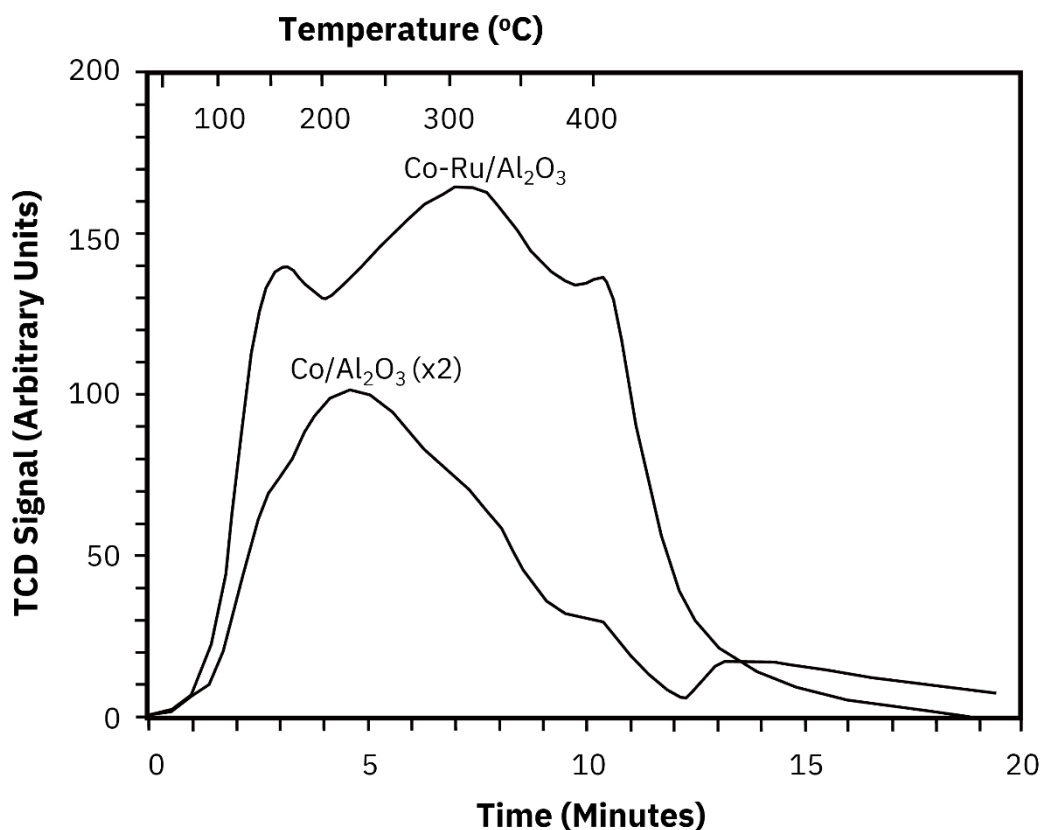


Figure 1: H_2 TPD curves of Al_2O_3 -supported Co catalysts, with and without Ru promoter; top x-axis shows desorption temperature, bottom x-axis shows heating time

provides valuable qualitative information about the strength of chemisorption based on the temperature distribution of the desorption profile.

Catalyst	Promoter	Method	H_2 Uptake ($\mu\text{mol/g cat.}$)
20 wt% Co/ Al_2O_3	None	Volumetric Chemisorption	41
20 wt% Co/ Al_2O_3	None	TPD	42
0.5 wt% Ru/20 wt% Co/ Al_2O_3	Ru	Volumetric Chemisorption	164
0.5 wt% Ru/20 wt% Co/ Al_2O_3	Ru	TPD	188

Table 2: Comparison of volumetric chemisorption and TPD H_2 uptake data for Al_2O_3 -supported Co catalysts with and without Ru promoter

4. Conclusions

This study demonstrates that both dynamic (pulse chemisorption and TPD) and static (volumetric) chemisorption techniques, when properly applied, yield consistent and reliable results across a variety of catalysts. Dynamic methods offer clear advantages in terms of simplicity, flexibility, and significantly reduced analysis time—often cutting hours down to minutes. Additionally, dynamic

techniques provide enhanced sensitivity and the potential for gaining qualitative insights into adsorption strength and surface interactions, making them ideal for both research and quality control environments.

The **AMI 300** and **AMI 400**

series of chemisorption analyzers provide highly precise and accurate measurements for static (volumetric) chemisorption, pulse chemisorption, and TPD experiments. They are also highly customizable, allowing for a variety of sample types, unique experiment design, and optional *in situ* spectrometry. These features, described in Figure 2, make **AMI** chemisorption analyzers an excellent tool for a broad range of chemisorption experiments.



Figure 2: Highlighted features of **AMI 300** and **AMI 400** series of chemisorption analyzers

5. References

- (1) Chien, A. C. and Chuang, S. S. C. Static and dynamic hydrogen adsorption on Pt/AC and MOF-5. *Int. J. Hydrogen Energy*, **2011**, *36*, 6022-6030.
- (2) Rioux, R. M.; Hoefelmeyer, J. D.; Grass, M.; Song, H.; Niesz, K.; Yang, P.; Somorjai, G. A. Adsorption and Co-adsorption of ethylene and carbon monoxide on silica-supported monodisperse Pt nanoparticles: Volumetric desorption and infrared spectroscopy studies. *Langmuir*, **2008**, *24*, 198-207.
- (3) Chantaravitoon, P.; Chavadej, S.; Schwank, J. Pt-Sn/Al₂O₃ catalysts: Effect of catalyst preparation and chemisorption methods on H₂ and O₂ uptake. *Chem. Eng. J.* **2004**, *98*, 99-104.
- (4) Reuel, R. C. and Bartholomew, C. H. Effects of support and dispersion on the CO hydrogenation activity/selectivity properties of cobalt. *J. Catal.* **1984**, *85*, 78-88.
- (5) Kogelbauer, A.; Goodwin, J. G. Jr.; Oukaci, R. Ruthenium promotion of Co/Al₂O₃ Fischer-Tropsch catalysts. *J. Catal.* **1996**, *160*, 125-133.